

CatchGene[®] FFPE Tissue RNA Kit

Cat. No.	Rxn		
MR21004	4		
MR21050	50		
MR21250	250		

Kit Description

The CatchGene[™] FFPE Tissue RNA kit provides a safer, quicker and more efficient silica membrane-based RNA purification from Formalin-fixed, Paraffin-embedded FFPE tissue sample. Traditional method in pre-treatment of FFPE tissue sample uses xylene, which is highly toxic and carcinogenic . Hence, CatchGene FFPE Tissue RNA Kit offers a safer and optimized xylene-free method for RNA purification from FFPE tissue. Furthermore, Buffer RFTL and RFB is able to lyse FFPE tissue efficiently, and eventually minimize vortex steps to prevent fragmentation and preserve high integrity of RNA.

Kit Storage

Kit Content

Kit Content						Kit Stolage	
		4rxn	50rxn	250rxn		Upon arrival,	
	MR21 Column	4	50	250	pcs	1. Please store MR21 Column at 4°C	
	Collection Tubes (2 ml)	12	150	750	pcs	for long term storage.	
	Buffer AE	0.5	1.5	10	ml	2. Please store Proteinase K at -20 $^{\circ}$ C	
	Proteinase K	1x2	11x2	11x10	mg	for long term storage.	
	Buffer DWX	2	27	135	ml	Buffer, solvent and consumables, please	
	Buffer RFTL	1.1	13.5	67.5	ml	store at 15-25 $^{\circ}$ C.	
	Buffer RFB	1.1	13.5	67.5	ml		
	Buffer RW1 (concentrated)	3.36	42	105x2	ml	If a precipitate has formed in Buffer RFTL	
	Buffer RW2 (concentrated)	0.68	8.4	42	ml	or RFB, dissolve by incubating at 60°C and	
	RNase-Free H ₂ O	0.96	12	60	ml	cool down to 25°C for using.	

Kit Preparation

- Prepare 20 mg/ml Proteinase K
 For 1 mg Proteinase K, please add 50 μl Buffer AE into tube and vortex thoroughly for dissolving.
 For 11 mg Proteinase K, please add 550 μl Buffer AE into tube and vortex thoroughly for dissolving.
 After dissolving into the solvent, please store at 4°C for 6 month or -20°C for 1 year.
- 3. Prepare Buffer RW1 Add equal volume of 100% EtOH into Buffer RW1 (concentrated) to get Buffer RW1. After adding 100% EtOH, please check the sticker on the bottle and close the cap tightly.
- Prepare Buffer RW2
 Add 4 volume of 100% EtOH into Buffer RW2 (concentrated) to get Buffer RW2.
 After adding 100% EtOH, please check the sticker on the bottle and close the cap tightly.

General Protocol

- 1. Place 5-10 μm sections (up to 4 sections) in the micro-centrifuge tube (not provided). Add 450 μl DWX buffer, vortex vigorously for 15 sec. Spin down to collect sample in the bottom.
- 2. Incubate at 60° C for 20 min. Brief spin down.
- 3. Add 225 μl Buffer RFTL (Please add 1% ß- mercaptoethanol freshly) and mix thoroughly by vortex 15 sec.
- 4. Centrifuge at 11,000 x g for 1 min.
- 5. Add 20 μl Proteinase K (20 mg/ml) to the lower clear phase. Mix gently by pipetting.
- 6. Incubate at 60° C for 30 min. (Or until all tissue lysed, longer incubation time will increase yield of some FFPE tissue.) Brief spin down.
- 7. Incubate at 80 $^\circ\!\mathrm{C}$ for 15 min.
- 8. Add 225 μI Buffer RFB in to the lower phase, mix gently by pipetting.
- 9. Centrifuge at 11,000 x g for 1 min.
- 10. Aspirate 400 μ l lower clear phase lysate into a new 1.5 ml micro-centrifuge tube.
- 11. Add 200 µl of 100% EtOH and mix thoroughly by vortex for 5 sec, brief spin down.
- 12. Transfer all mixture to Spin Column (with 2ml Tube), Centrifuge at 11,000 x g for 1 min. Discard the flowthrough and change a new Collection Tube.
- 13. Add 700 µl Buffer RW1 into spin column, centrifuge at 11,000 x g for 1 min, discard the flow-through.
- 14. (Optional) On column digest of DNA with DNase I (not provided).
- 15. Add 700 μl Buffer RW1 into spin column, centrifuge at 11,000 x g for 1 min, discard the flow-through.
- 16. Add 700 μl Buffer RW2 into spin column, centrifuge at 11,000 x g for 1 min, discard the flow-through.
- 17. Change a new Collection Tube, centrifuge at 11,000 x g for 3 min.
- 18. Place the spin column into 1.5 ml Elution Tube, add 30-100 μl RNase-Free H_2O and incubate at 25°C (room temperature) for 3 min.
- 19. Centrifuge at 11,000 x g for 1 min for elution.

FOR RESEARCH USE ONLY